

# RAPID EPS

## RAPID EPS Instruction Manual

Thank you for purchasing our RAPID EPS.  
Please read this manual carefully before use and  
save this manual for your future use.

Please note that product specifications and  
packings may change without prior notice.

**Announcement for your use with ease.**

### 【Purpose of use】

This product can be used for research purposes only.

### 【Notes for use】

Handle this product carefully when unpacking from the bag.

Do not process this product in autoclave treatment.


### 【Storage and disposal of product】

Store at room temperature and in a place not exposed to direct  
sunlight. To dispose the product after use, comply with local laws  
and ordinances, and process properly according to the regulations  
of your facility.

## Instructions


\*Please wear gloves to handle the product.

1

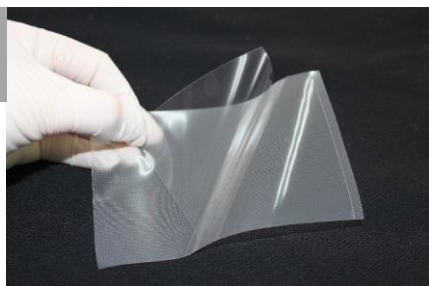

Remove the release liner from the slit along the edge  
and expose the adhesive surface.

2

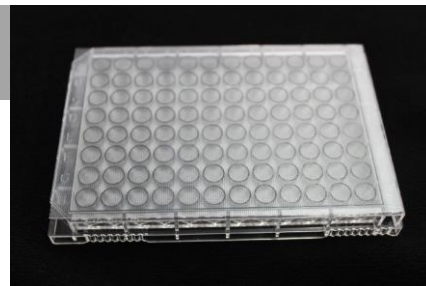

Adhere the adhesive surface on a well plate.  
After pasting, press the surface with hands or a  
spatula, etc. and **adhere it firmly.**

## Manufacturer



BioChromato, Inc.

1-12-9 Honcho, Fujisawa, Kanagawa-ken 251-0053 JAPAN

TEL : +81-466-23-8382 FAX : +81-466-23-8279

E-mail : [oversea@bicr.co.jp](mailto:oversea@bicr.co.jp)

Website: <https://biochromato.com/>

Catalog Code	REPS001 R37.001.00
Product Name	RAPID EPS
Bonding method	Pressure sensitive adhesive
Materials	polyolefin
Operating temperature limit	-80°C~80°C
Size	W 80 mm × L 122 mm
Quantity	100sheets/Box